

Reports the Way You Need Them from a Searchable Database

- Custom Report Generation from ScintiVision-32
- Multiple custom report formats via specific templates
- Spectral plot included with report
- Creates Access®-compatible results database
- Archival capability

**For Gamma Spectroscopy
with NaI Detectors**

Reports Your Way Report Writer (A46-B32), in conjunction with ScintiVision-32 software for NaI detectors, lets you produce a wide range of output reports — in exactly the form you require.

Instantly Searchable Database Report Writer reads in the analysis results from the ScintiVision-32 output file, stores them into an Access-compatible database, and produces the report. Thereby, you progressively generate a PC-searchable database of all your analysis results. Then use Access to quickly generate, for example, daily summaries, exception reports, and cumulative release reports. Report Writer thus acts as a bridge between ScintiVision-32 results and database applications or a laboratory information management system.

Our Template or Yours The output report is generated using a template file. The default template file supplied reproduces the standard ScintiVision-32 output report. Additional user-defined files¹ may be used to produce totally custom output reports via this template development tool.

In addition to automatic output to the printer, the reports may be saved to disk in various industry-standard formats:

ASCII, Rich Text, Crystal Reports, WORD, Comma-separated.

Multicolor Plots Automatically A plotting program is included which permits the spectrum to be plotted to the printer automatically as part of the output report. The plot attributes may be prespecified.

As Automatic as You Want to Be Report Writer can be run independently in order to regenerate or reformat reports from previous analyses, or automatically as part of the ScintiVision-32 analysis sequence, in which case the standard internal ScintiVision-32 output report is disabled.

Good Housekeeping A facility is provided to allow database and supporting files to be archived to a specified directory. File types may be chosen from templates, spectral files, output files, and region-of-interest files. The archival process requires only a single click. Similarly a “purge” option allows removal of such files to the system recycle bin or permanent deletion.

¹In Crystal Report Writer, V5.0 format; trademark of Crystal Software, Vancouver, BC.

Report Writer for ScintiVision™-32 A46-B32

Ordering Information

To order, specify:

Model	Description
A46-B32	Report Writer option for ScintiVision-32
A46-U32	Update to A46-B32
A46-G32	Documentation for A46-B32

Prerequisites

A46-B32 will function satisfactorily on any system capable of supporting ScintiVision-32

For custom report template development, a copy of Crystal Reports Version 5.0 or later is required.

Specifications subject to change
011007

ORTEC®

www.ortec-online.com

Tel. (865) 482-4411 • Fax (865) 483-0396 • info@ortec-online.com
801 South Illinois Ave., Oak Ridge, TN 37831-0895 U.S.A.
For International Office Locations, Visit Our Website

AMETEK®
ADVANCED MEASUREMENT
TECHNOLOGY