



ORTEC®
Sunpower
Princeton Applied Research®
Solartron Analytical®
Signal Recovery®



ADVANCED MEASUREMENT TECHNOLOGY
801 South Illinois Avenue, Oak Ridge, TN 37830
Tel: 865.482.4411 Fax: 865.483-0396

GAMMAVISION PRODUCT REGISTRATION

Steps to generate the ORTEC Registration Form

GammaVision requires product registration within 60 days after installation. Until registered, the dialog shown below opens each time the program is started. Once the Registration key is obtained, it can be entered in the "Registration Key" box within the dialog shown below and confirm registration by clicking the "Register Now" button. Once this is complete, the dialog will no longer be displayed and the software will be properly registered.

GammaVision Registration

Installation Key: 056-5885534

Registration Key:

Registration Information

Please register the software by e-mail at ORTEC.Software@Ametek.com, or call ORTEC at extension #449 to obtain a registration key. When you register, please have your Registration Form ready.

Days Left: 60

Generate ORTEC Registration Form

Register Later Register Now

To complete the registration form and submit to ORTEC to obtain the required registration key, please complete the following steps:

1. Click Generate ORTEC Registration Form to open the ORTEC Registration Form dialog box.
2. Enter the Serial Number for the copy of GammaVision purchased. The serial number is located on the box.
3. Click the Write ORTEC Registration File button to create a text file containing the registration information. A standard file-save dialog will open. Enter a filename and location, then click SAVE. Email this file to the ORTEC email address listed on the GammaVision Registration dialog. ORTEC will respond with the registration key.
4. Enter the registration key on the GammaVision Registration dialog and click Register Now.